

DS2500 Solid or Liquid Analyzer

Preventative Maintenance Guidelines



RECOMMENDED MAINTENANCE

Regular maintenance at defined intervals is required to keep your DS2500 Solid or Liquid Analyzer operating at peak performance so that your results stay accurate and trustworthy, and you minimize instrument downtime.

Metrohm USA recommends several simple actions to ensure your instrument is always ready for operation. Our recommended schedule is below:

Action	Note*
Fan filter replacement	Monthly, or as needed
Lamp exchange	3-6 months, or as needed
Cooling fluid refill	Annually, or as needed
Cleaning of exterior surfaces	Monthly, or as needed
Cleaning of glass surfaces	As needed with brush or microfiber cleaning cloth when dirty

* Maintenance schedule is dependent on environmental conditions and usage

RECOMMENDED CONSUMABLES AND SPARE PARTS

Consumable and spare parts can be ordered online at usa-shop.metrohm.com or by calling Metrohm USA customer support at 866-METROHM.

Note: Have isopropanol and optical wipes available to clean surfaces and glass parts of the analyzer.

Product	Part Number
DS2500 fan filter - 5 pack	49220490
DS2500 single fan filter with holder	49220300
DS2500 Spare Lamp	67430050
DS2500 Cooling Fluid	49220530
Microfiber cleaning cloth	67430060
Brush	SNG-NIR036778

**ADDITIONAL RECOMMENDATIONS FROM YOUR
FIELD SERVICE ENGINEER:**

ACCESSING TECHNICAL SUPPORT

Technical support is available by emailing techsupport@metrohmusa.com or calling 1-866-METROHM from 8:30AM to 7PM EST, Monday-Friday.