

MagIC Net – Exporting Determination Files

MagIC Net – How to Export .idet files


Technology: Ion Chromatography

Overview


There are 2 steps necessary to export data files. First, an export template must be created. Once the template is created, the .idet data files can easily be exported from the software.

Instructions for Template Creation

1. Go to Database > Tools > Templates > Select Export Templates

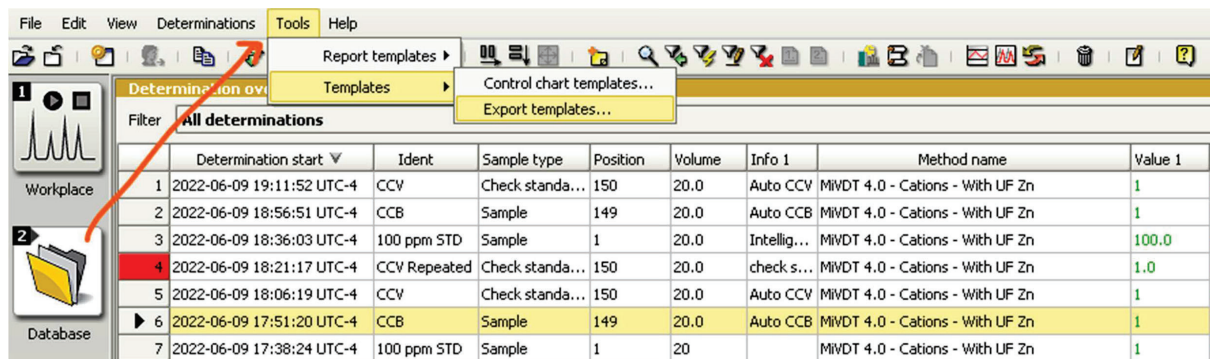


2. The Export Templates Window opens. Select "New".
3. The Export Template "New File" Window will open.
4. Rename and create a comment if desired.
5. Select the target directory you wish to have the files exported to by pressing the
6. button. E.g., Flash drive, Desktop folder.
7. 6. File type = *.idet (MagIC Net format)
8. 7. File Name = Sample identification, Ident

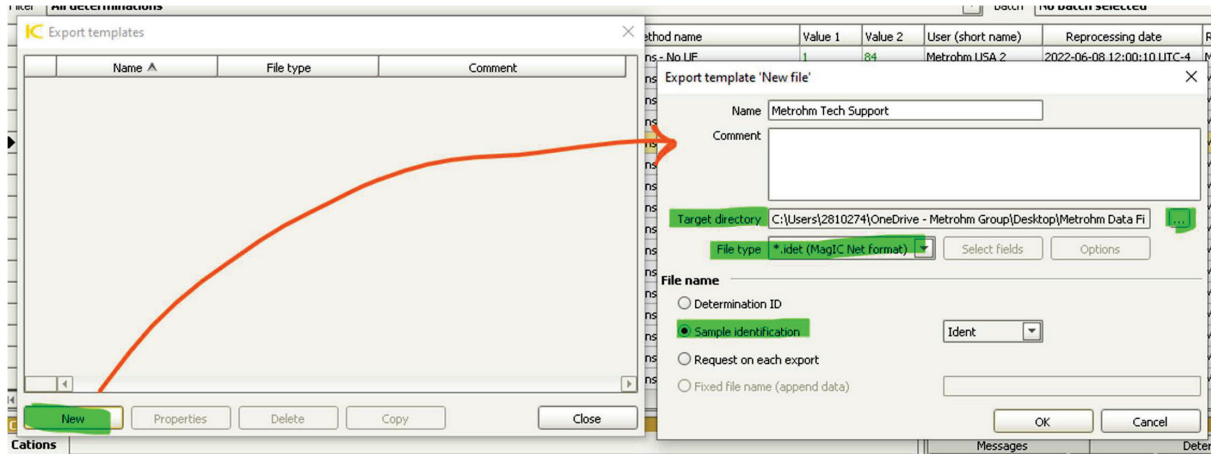


8. Select Okay to close out of the Export Template 'New File' window.

9. Select Close out of the Export Templates.

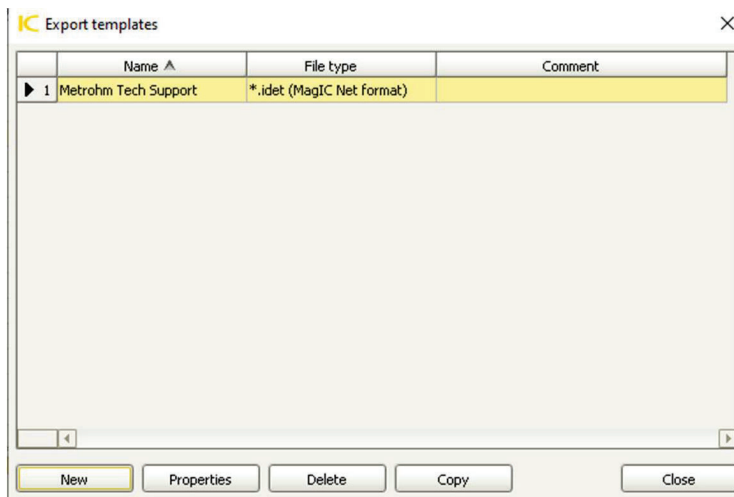


Instructions for Exporting Data Files


1. Go back into the database. Select the files you wish to export - by using CRL + click to
2. highlight all the files.
3. Select Determinations in the Upper Ribbon and choose Export...

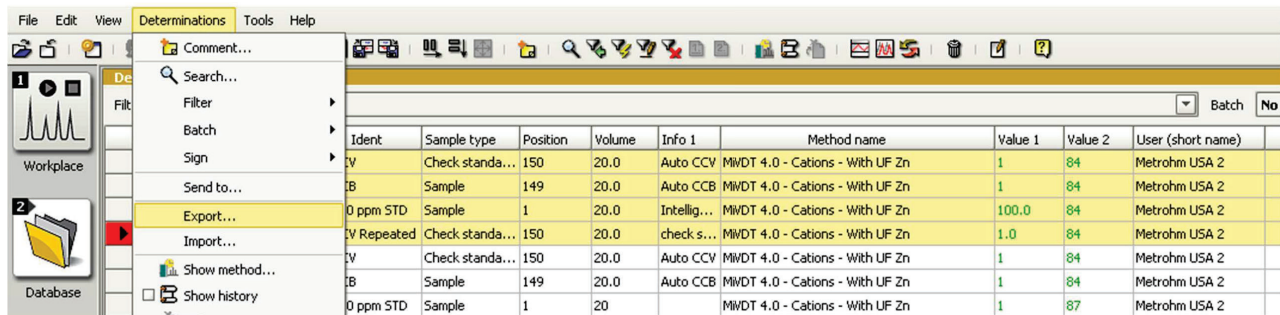


- 'Export Determinations' window will appear. Selection = All selected data records. Use the drop-down menu to select the Export Template previously created if it has not already been populated. Select 'OK'.

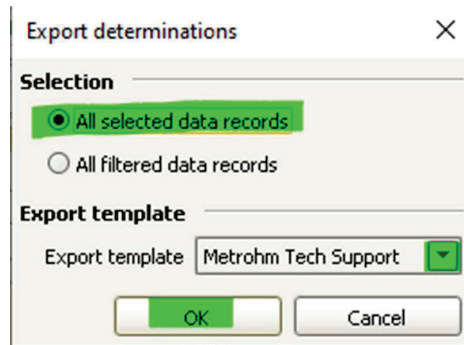

- The lower right-hand corner will have a status bar indicating the export is in progress.

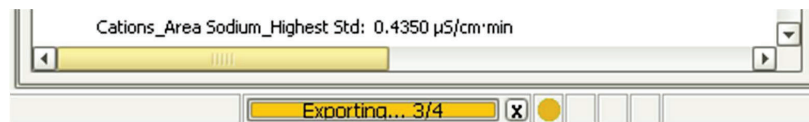

- Once complete, navigate to the folder designated for creating the export template to view the data files.

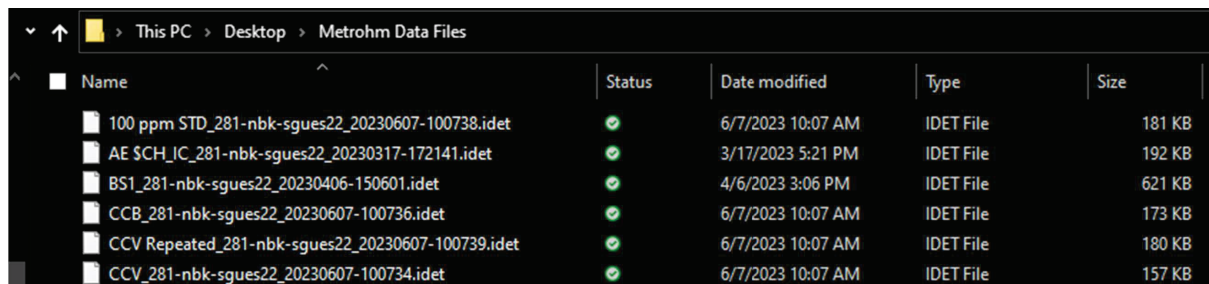

Name	Status	Date modified	Type	Size
100 ppm STD_281-nbk-sgues22_20230607-100738.idet	✓	6/7/2023 10:07 AM	IDET File	181 KB
AE \$CH_IC_281-nbk-sgues22_20230317-172141.idet	✓	3/17/2023 5:21 PM	IDET File	192 KB
BS1_281-nbk-sgues22_20230406-150601.idet	✓	4/6/2023 3:06 PM	IDET File	621 KB
CCB_281-nbk-sgues22_20230607-100736.idet	✓	6/7/2023 10:07 AM	IDET File	173 KB
CCV Repeated_281-nbk-sgues22_20230607-100739.idet	✓	6/7/2023 10:07 AM	IDET File	180 KB
CCV_281-nbk-sgues22_20230607-100734.idet	✓	6/7/2023 10:07 AM	IDET File	157 KB

- If necessary, move the exported files to a USB flash drive if email access is not on the MagIC Net PC.
- Zip the files up in a folder if sending to Metrohm Tech Support for review.